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**WWW.COMPRABASCULA.COM**

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## 1. INTRODUCTION

**NOTE:** These balances are not apt for the uses indicated in the section second of the directive 90/384/CE.

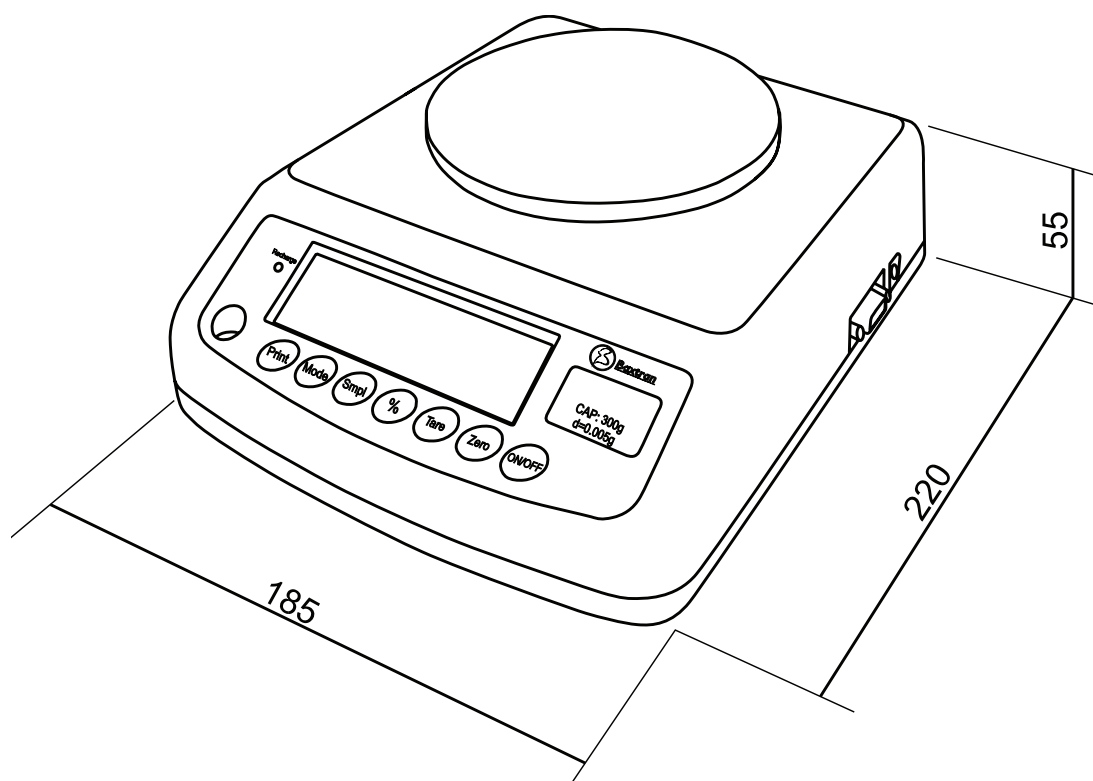
The BAT series of electronic balances provides an accurate, fast and versatile series of general purpose balances with counting and % weighing functions.

There are 4 models in this serie, with capacities up to 3000g.

They all have stainless steel weighing platforms on an ABS base assembly.

All the keypads are light touch switches and the displays are large easy to read liquid crystal type displays (LCD). The LCD's are supplied with a backlight and have loading parts counting.

All units include automatic zero tracking, automatic tare, and an accumulation facility that allows the count to be stored and recalled as an accumulated total.

## 2. SPECIFICATIONS

### 2.1 SERIE BAT

SERIE BAT				
N° Model	BAT300	BAT600	BAT1500	BAT3000
Capacity	300	600	1500	3000
Readability (g)	0,005	0,010	0,020	0,050
Resolution	1/60.000	1/60.000	1/60.000	1/60.000
Tare range (g)	100	300	600	1000
Minimum capacity (g)	0,1	0,2	0,4	1,0
Linearity ± (g)	0,01	0,02	0,04	0,10
Pan size	Ø 120 mm		136 x 162 mm	
Wind protector	SI	SI	NO	NO
Units of measurement	g / ct / lb / oz / d / gn / ozt / dwt / mm / tl. T / tl. c / tl. t / t.			

### 2.2 COMMON SPECIFICATIONS

Output	Optional RS.232 output
Stabilization time	Habitually 2 seconds
Operating temperature	0 °C - 40 °C / 32 °F - 104 °F
Power supply (external)	9V / 800 mA o rechargeable internal battery
Calibration	Automatic external
ACDC	$\Sigma$ - $\Delta$
Internal divisions	1.000.000
Screen	LCD 6 digits; 20 mm with backlight and bar of charge.
Materials	ABS, stainless steel pan
Pan size	Ø 120 mm / 136 x 162 mm
General dimensions (wide x width x height)	180 x 220 x 85 mm
Gross weight	1,4 kg (it includes the rechargeable battery)
Other components and specifications	High accuracy in the counting way and percentage weight

## 3. INSTALLATION

### 3.1. GENERAL INSTALLATION

The scale should be sited in a location that will not degrade the accuracy.

Avoid extremes of temperature. Do not place in a direct sunlight or near air conditioning vents.

Avoid unsuitable tables. The tables or floor must be rigid and not vibrate. Do not place near vibrating machinery.

Avoid unstable power sources. Do not use near large users of electricity such as welding equipment or large motors.

Avoid air movement such as from fans or opening doors. Do not place near open windows.

Keep the scale clean.

This scale has not been designed like a waterproof scale (Ip44) reason why must avoid high humidity levels that could cause condensation. Avoid the direct bonding with the water. Do not sprinkle water on the scale and do not submerge it. If the scale makes contact with water, the reading data can be unstable, or the scale can not work correctly, in that case, extinguishes the electrical power immediately.

Do not stack material on the scales when they are not in use.

### 3.2. INSTALLATION OF BAT SERIES

Level the scale by adjusting the four feet. The scale should be adjusted such that the bubble in the spirit level is in the center of the level and the scale is supported by all four feet. If the scale rocks readjust the feet.

Attach the adapter to the connector on the side of the scale. Please use the power as the label.

Mount the protective bell with the five plastic pieces included in all the models except those of 1.5 kg and 3 kg of capacity.

Place the protective bell so that is fitted perfectly on the scale and on does not move, downwards leaving small notches located at corners of the bell.

Do a calibration of weight as it is described at section 8 (Configuration) in the function Normal Calibration (P2CAL).

## 4. KEYPAD DESCRIPTION

KEY	DESCRIPTION	
	PRIMARY FUNCTION	SECONDARY FUNCTION
<b>Zero</b>	Set the zero point for all subsequent weighing. The display shows zero.	"Enter" key when setting parameters or other functions.
<b>Tare</b>	Tares the scale. Stores the current weight in memory as a tare value, subtracts the tare value from the weight and shows the results. This is the net weight.	
<b>%</b>	Enters the percent weighing function. When in percentage mode and current value is zero, press this key to return normal weighing mode.	
<b>Print</b>	To print the results to a PC or printer using the optional RS-232 interface. It also adds the value to the accumulation memory if the accumulation function is not automatic.	"ESC" is to return to normal operation when the scale is in a parameter setting mode.
<b>Mode</b>	This key will select either kilograms, pounds, ounce for the weighing unit when weighing mode. When counting mode, allows the weight, unit weight, and count to be seen when parts counting.	A secondary function, change current value for parameters or other functions.
<b>Smpl</b>	Use this key to enter counting mode when normal weighing mode.	Use this key to return normal weighing mode when counting mode.
<b>On/Off</b>	Use this key to turn on or turn off the power.	

## 5. OPERATION

### 5.1. ZEROING THE DISPLAY

You can press the **Zero** key at any time to set the zero point from which all other weighing and counting is measured, within 10% of power up zero. This will usually only be necessary when the platform is empty. When the zero point is obtained the display will show the indicator for zero.

The scale has an automatic rezeroing function to account for minor drifting or accumulation of material on the platform. When the weight goes over 20 % of the full capacity, "read zero" function is obsolete.

### 5.2. TARING

Zero the scale by pressing the **Zero** key if necessary. The zero indicator will be on.

Place a container on the platform, a value for its weight will be displayed.

Press the **Tare** key to tare the scale. The weight that was displayed is stored as the tare value and that value is subtracted from the display, leaving zero on the display. The "TARE" indicator will be on. As product is added only the weight of the product will be shown. The scale could be tared a second time if another type of product was to be added to the first one. Again only the weight that is added after taring will be displayed.

When the container is removed a negative value will be shown. If the scale was tared just before removing the container this value is the gross weight of the container plus all product that was removed. The zero indicator will also be on because the platform is back to the same condition it was when the **Zero** key was last pressed..

### 5.3. WEIGHING A SAMPLE

To determine the weight of a sample first tare the empty container then place the sample in the container. The display will show the weight and the units of weight currently in use.

### 5.4. PERCENT WEIGHING

The scale will allow a sample weight to be shown as 100%. Then any other weight placed on the scale will be displayed as a percentage of the original sample. For example is 350g is placed on the scale and the **%** key is pressed the display will show 100.00%.

Removing the 350g weight and putting a 300g weight on the scale the display will show 85.71% as 300g is 85.71% of 350g.

Note: the scale may jump by large numbers unexpectedly if small weights are used to set the 100% level. For example if only 23.5g is on a scale with 0.5g increments and the scale is set to 100%, the display will show 100.00%, however a small change of weight will cause the display to jump to 102.13% as one scale division (0.5g) increase to 24.0g will be equivalent to a 2.13% increase.

Removing the 350g weight and putting a 300g weight on the scale the display will show 85.71% as 300g is 85.71% of 350g.

Pressing the **%** key will return the scale to weighing.

## 5.5. PARTS COUNTING

Before beginning, tare the weight of any container that will be used, leaving the empty container on the scale. Place the number of samples on the scale. The number should match the options for parts counting, 10, 20, 50, 100 or 200 pieces.

When the scale is showing weight, pressing the **Smpl** key will start the parts counting function.

The scale will show "SP 10" asking for a sample size of 10 parts. Pressing the **Mode** key the display will cycle through the options: 10, 20, 50, 100, 200 and back to 10.

Press the **Smpl** key again when the number matches the number of parts used for the sample. As more weight is added the display will show the number of parts (pcs).

Press the **Mode** key to display unit weight (g/pcs), Total weight (kg) or the count (pcs).

Press the **Smpl** key to return to normal weighing.

## 5.6. ACCUMULATION

When the scale is set to manual accumulation the weight displayed will be stored in memory when the **Print** key is pressed and the weight is stable.

The display will show "ACC 1" and then the total in memory for 2 seconds before returning to normal. If the optional RS-232 interface is installed the weight will be output to a printer or PC.

Remove the weight, allowing the scale to return to zero and put a second weight on. Press **Print**, the display will show "ACC 2" and then show the new total. Continue until all weights have been added.

To view the totals in memory press the **Print** key when the scale is at zero. The display will show the total number of items "ACC xx" and the total weight before returning to zero. The totals will also be printed via the RS-232 interface.

To erase the memory press **Print** to view the totals and then press the **Smpl** key to clear the memory.

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When the battery needs to be recharged, the indicator  lights. Please, load the battery or use an adapter of alternating current when he/she sees this lit indicator.

The scales can work with batteries if one wants. The duration of the battery is of approximately 72 hours.

To load the battery plug it to the electric net. It is not necessary to light the scale.

The recharge should last 12 hours to fill the battery to full capacity.

There is a LED that indicates the state of the load of the battery. When the scale is plugged to the electric net, the battery goes into it will load. If LED is green the battery has complete load. If it is red, the battery needs load, and the yellow indicates that the battery is in recharge process.

With the use, the battery can go losing capacity to make complete loads. If the duration of the load of the battery becomes unacceptable, contact its distributor.

Note: the new batteries are partially loaded. Before being able to use their scale, it should install and to load the battery following the following instructions.

Some batteries have a better yield after several complete cycles of load / it discharges.

The yield of the batteries depends on numerous factors, included the configuration of the retroiluminación and the operative one.

Not never use a loader or a battery that are damaged.

Not make a short circuit in the battery. A short-circuit can take place accidental when an object metallic (currency, clip or pen) cause one direct connection of the poles + and. of the battery (you throw metallic of the battery), for example when it takes a reserve battery in the pocket.

To make a short-circuit of the poles can damage the battery or the object that it is connected.

Not throw the batteries to the fire.

Throw the batteries following the local (for example, reciclaje) regulation.

Not throw the batteries like domestic garbage.

Avoid to load the battery under conditions of lack of air.

In order to maximize the yield of their battery:

Always use batteries and adapter of CA originals. The guarantee of the scale doesn't cover the damages caused as consequence of the use of other batteries and/or different loaders.

The tension of exit of adapting CA is of 9V, but the range of normal tension will oscillate between the 11V and the 15V.

The new batteries or those batteries that have been stored during long periods of time can require a bigger time of load.

Maintain the battery to ambient temperature or a next temperature when it makes the load.

Not expose the batteries to inferior temperatures to -10°C or superiors at 45°C.

In the course of periods of long time, the batteries go losing load capacity progressively and they require longer times of recharge.

This is normal. If you load the battery regularly and he/she observes that it diminishes the period of operation or that it increases the period of load, it is probably moment of acquiring a new Battery.

## 7. RS.232 OUPUT

The BAT series scales can be ordered with an optional RS-232 output.

Specifications:

- RS-232 output of weighing data
- ASCII code
- 9600 Baud
- 8 data bits
- No Parity

Connector: 9 pin d-subminiature socket

- Pin 3: Tx (Output)
- Pin 2: Rx (Input) (not used at this time)
- Pin 5: GND

BAT scales can be connected with PC, remote display or mini printer, setup see section 8.

## 8. PARAMETERS

The scale has 8 parameters that can be set by the user plus a method of entering the calibration section. To set parameters press the **Mode** key during self checking.

The display will show the first function, "F1 UNIT". Pressing the **Mode** key will cycle through the other functions.

Pressing **Zero** key will allow you to set the function. Use the **Print** key to leave a parameter unchanged.

Note: when display shows "TECH", press **Zero** key to enter, display shows "PIN", press **Smpl** , **Print** , **Tare** key, and press **Zero** key to enter technical parameters setting mode.

## PARAMETERS TABLE

FUNCTION	DESCRIPTION
F1 UNT	Sets the which units want to use: g / ct / lb / oz / d / gn / ozt / dwt / mm / tl.T / tl.c / tl.t / t use <b>Mode</b> to switch, use <b>Zero</b> key to sure.
F2 EL	Set the backlight to be on, automatic or off. EL on: Backlight always on EL Au: Backlight automatically turn on EL off: Backlight allways off Use <b>Mode</b> to switch, use <b>Zero</b> key to sure.
F3 PRT	P cont, continuous sending. P auto, when the weight is stable sends the data. P prt, printer. Press <b>Zero</b> , now you can select the speed of transmission (600, 1200, 2400, 4800 or 9600 bps), press <b>Mode</b> to change speed and <b>Zero</b> to save the wished speed. Now it selects the printer (LP-50 or TP), press <b>Mode</b> to change printer and <b>Zero</b> to save.
TECH	Pin Enter the programming and calibration menus by entering the correct password ( <b>Smpl</b> <b>Print</b> <b>Tare</b> ) and press <b>Zero</b> key to enter.
P1 LIN	Linear calibration. Press <b>Zero</b> . Display will show PIN again. Press keys <b>Print</b> , <b>Smpl</b> , <b>%</b> y <b>Zero</b> to enter. Display will show UNLOAD. Take off all the weight of the tray. Display will show the first request of weight for the calibration. Place this weight on the tray, when the balance is stable ignites the stability indicator and accepts this weight automatically. After the first calibration of weight, display will show the following calibration weight. Place this weight on the tray, when the balance is stable ignites the stability indicator and accepts this weight automatically. After the second calibration of weight, display will show the following calibration weight. Place this weight on the tray, when the balance is stable ignites the stability indicator and accepts this weight automatically. The scale will begin to count backwards until reaching zero after it has accepted the last weight. During the count down, you must retire all the weight of the platform. The scale becomes in normal weighing mode.
P2 CAL	After entering in the calibration mode, display will show UNLOAD. Take off all the weight of the tray. When the scale is stable, it will fix initial zero point automatically. Then display will show LOAD. Place the weight for the calibration on the scale. Minimum calibration weights: BAT300, 50g. BAT600, 100g. BAT1500, 300 g. BAT3000, 500 g. Place this weight on the tray, when the balance is stable ignites the stability indicator and accepts this weight automatically. The scale will begin to count backwards until reaching zero after it has accepted the weight. During the count down, you must retire all the weight of the platform. The scale becomes in normal weighing mode.
P5 GRA	Gravity. It appears the value of gravity, to change this value use the keys: <b>Mode</b> , places the value to 0. <b>Tare</b> , increases the value of the active digit. <b>Smpl</b> , changes the active digit at right. <b>%</b> , changes the active digit at left. Press <b>Zero</b> to save.

## 9. APPENDIX

### 9.1. ERROR CODES

During the initial power-on testing or during operation it is possible the scale may show an error message. The meaning of the error messages is described below.

If an error message is shown repeat the procedure that caused the message, turning the balance on, calibration or other functions. If the error message still is shown then contact your dealer for further support.

ERROR CODE	DESCRIPTION	POSSIBLES CAUSES
<b>ERR 4</b>	Initial Zero is greater than allowed (4% of maximum capacity) when power is turned on or when the <b>Zero</b> key is pressed.	Weight on the pan when turning the scale on. Excessive weight on the pan when zeroing the scale. Improper calibration of the scale. Damaged load cell. Damaged electronics.
<b>ERR 5</b>	Keyboard error.	Improper operation of the scale.
<b>ERR 6</b>	A/D count is not correct when turning the scale on.	Platform not installed. Load cell damaged. Electronics damaged.

To see the divisors from A/D converter, push **Zero** key and **%** key at once while the scale is doing the initial revision when it is turning on. Push **Zero** key to return at normal function.

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Declare under our responsibility that the denominated product:

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To which declaration refers to, conform with the followings standards or other normative documents:

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20049/108/CE directive on the electromagnetic compatibility.

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And the following harmonized norms/ normes harmonisées:

EN55022: 2006+A1:2007

EN61000-4-2, 4-3, 4-4, 4-5, 4-6, 4-11

EN55024:1998+A1: 2001+A2: 2003

EN61558-2-6:1997

Jordi ribalta

Director General  
General Manager

***Baxtran***